NOTICE TURBOTEST 33/300 AA


FEUILLE/SHEET 1/19

Code: 51006921

Machine: 1996-002

230V-50Hz

Révision du 25/09/1996

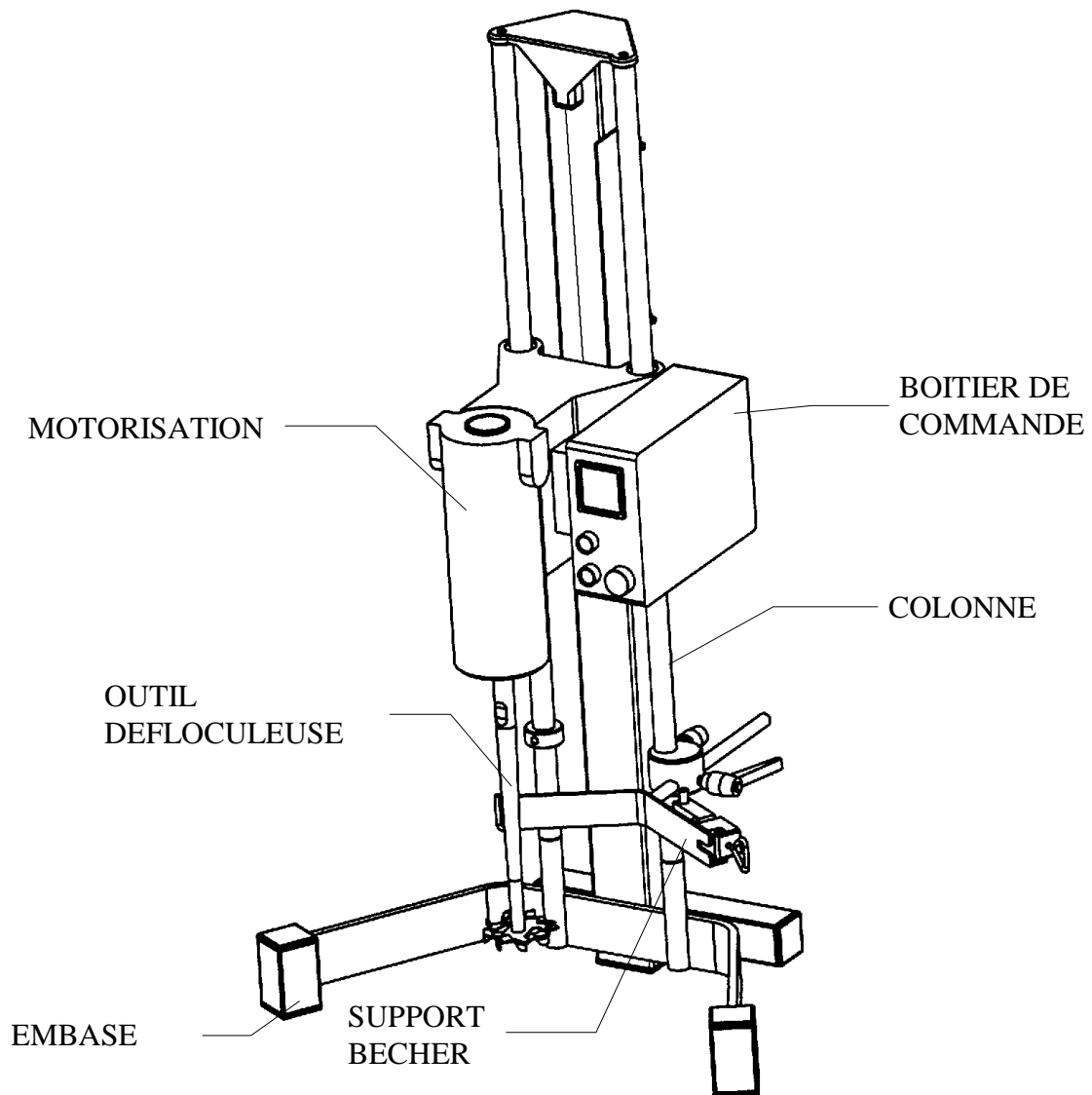*FIGURE NON CONTRACTUELLE
PICTURE NOT CONTRACTUAL*

Boitier de commande	Control panel
Colonne	Column
Embase	Base
Motorisation	Motorization
Outil défloculeuse	Defloculating propeller tool
Support becher	Beaker stay

NOTICE TURBOTEST 33/300 AA

Code: 51006921

Machine: 1996-002

230V-50Hz

FEUILLE/SHEET 2/19

Révision du 25/09/1996

SOMMAIRE

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SUMMARY

1. GENERALITIES	3
1.1- Properties	3
1.2- General features	3
1.3- Size.....	4
1.4- Construction.....	5
1.5- Warnings	5
2. IMPLANTATION	5
2.1- Prior arrangements	5
2.2- Packing and Handling.....	6
3. COMMISSIONING	6
3.1- Setting up.....	6
3.2- Control panel.....	6
3.3- How to operate.....	7
3.4- Emulsifier assembling	7
3.5-The turbines and their apply	8/9
4. ADVICE FOR UTILIZATION	10
4.1- Safety instructions.....	10
4.2- Contra-indication	10
5. MAINTENANCE	10
6. TROUBLE SHOOTING	11
7.MECHANICAL DRAWINGS	12
7.1- General drawing	12
7.2- Locking detection drawing	13
7.3- Slide drawing	13
7.4- Pulley drawing	14
7.5- Safety lifting drawing	14
7.6- Nomenclature	15
8. ELECTRIC DRAWINGS	17
8.1- Power supply 230V	17
8.2- Nomenclature	18
9. ANNEXE	19
9.1- Conformity	19
9.2- Extract from the conditions of sale	19
9.3- After sales service	19

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1. GENERALITIES

1.1- Properties

The TURBOTEST is a mixer designed to be used in laboratory. Its varied usage, so on realized mixes that on size of Beaker used, as well as the addition of products during the mix, necessitates its job by the personnel qualified. Its robust and simple idea, and its ergonomic stay to power lifting, make some an easy machine to manipulate. The large variety of turbines (more than 15 patterns) fact of the TURBOTEST a universal mixer for formulations in laboratory. Its silent functioning and its control panel render it usable in white hall. VMI-RAYNERI is not responsible for other utilizations or physical changes, except if we give prior written agreement.

1.2- General features

- Turbotest unloaded weight: 29,5 kg (with spindle + defloculating propeller turbine)
- Tool speed: 300 to 3300 rpm
- Motor power: 300 W
- Power supply: 230 V
- Frequency: 50 Hz
- Supply one phase + ground
- Noise level: 65 dBA max
- Protection of the control panel: IP54
- Beaker strap securing with beaker presence detector
- Total stroke: 380 mm
- Spindle diameter: 18 mm
- Spindle length: 260 mm
- Beakers size:
Diameter min: 80 mm / max: 220 mm
Height min: 100 mm / max: 300 mm
- Mixing capacity:
Min: 0,5 litres
Max: 10 litres

Nota: These capacities are given to indicative title and can vary following characteristics of products (viscosity, density...).

- Available accessories on order:
Emulsifier (Rotor-Stator type)
Stainless steel stay for 5 turbines
Electronic potentiometer with numeral speed display

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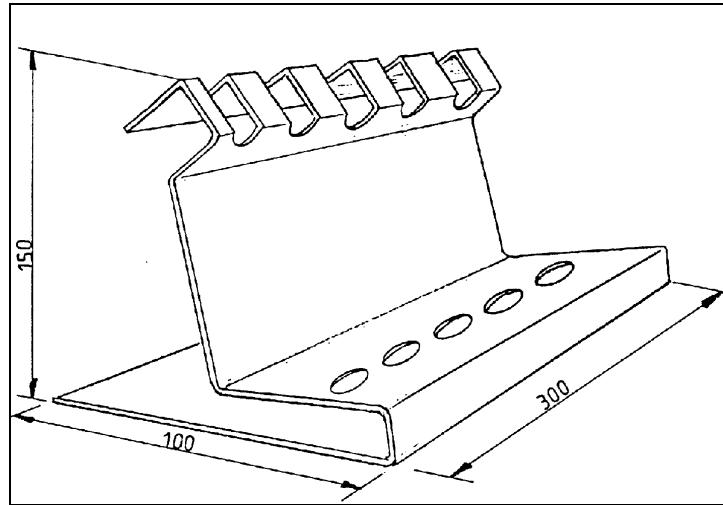
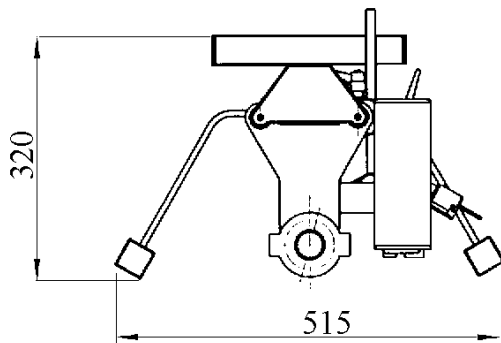
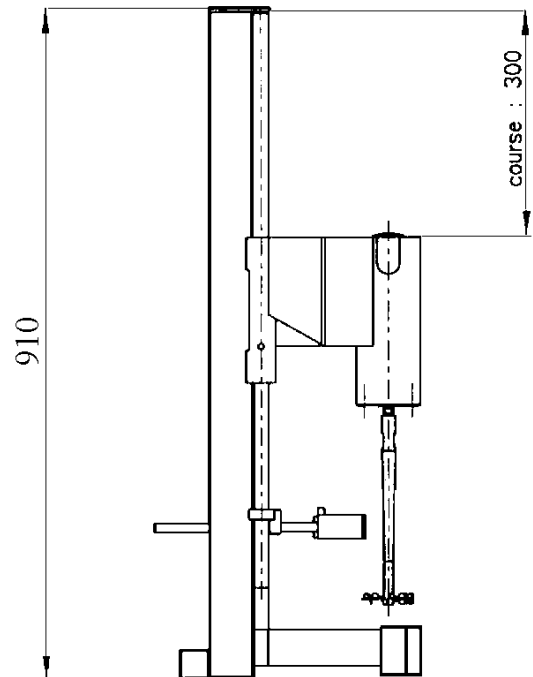
1.4- Construction

- Stay in stainless steel
- Columns of year guidance steel chromium-plated
- Motor slide in coated aluminium alloy of a paint polyurethane white RAL 9010
- Engine to continuous current and permanent magnets
- Spindle of agitation, fixing by cones and thread, in stainless steel AISI 316L
- Turbine in stainless steel AISI 316L polished

1.5- Warnings

- Every maintenance, repair or overhauling operation must be performed with duly qualified people with the machine disconnected from the power supply.
- It is strictly forbidden to introduce in the bowl all type of object useless to the product preparation.
- This operating manual must be kept carefully in order to make required maintenance. For any further request, remind the machine serial number (see identity plate fixed on the mixer side).

Information given within this manual may change and improve without any notice.

2. IMPLANTATION

2.1- Prior arrangements

- Inland packing:
L= 1000mm; l= 600mm; H= 350mm, W= 40 kg
- This machine is designed to be installed on a pallet, or any other plan to work.
- Release a surface on a length of approximately 800 mm and on a depth of approximately 600 mm minimum.
- Watch to a lighting sufficing the future posts work.
- The plan of laying has to be perfectly plan and horizontal, solid, stable and not to generate vibrations during the placement on the way.
- In the other case, for example a pallet tiled, it will be necessary to stall the support stable and durable.
- If the turbotest is not installed them the delivery, it will have to be stocked in its wrapping of origin, in a sheltered and dry place and to a room temperature.
- To anticipate a plug of feeding current, comprising a plug of earth adapted to the connection of masses, within 1 metre of the machine.

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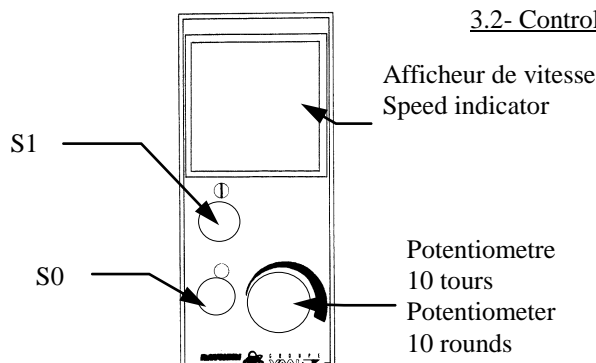
3. COMMISSIONING

3.1- Setting up

- Check that the power supply of the turbotest (230 V-one phase) indicated on the back face of the panel corresponds to the network supply.
- Unscrew the Bakelite plug placed on the left side on the slide, and lift the slide in high location.
- Lift the spindle with the 2 end wrenches supplied (key of 10 and 14).
- Adjust the position of the fixed Beaker with the 2 bakelite handles.
- Connect the turbotest to the power supply with the cable supplied (length 2 meters).
- Supply cable with the plug (16 A), phase, wire and ground.

Nota:

For special using, you can adjust the turbine height in low position by adjusting the stop placed on the left column, with the 6 sides wrench supplied in the machine.

For the turbine dismantling, use the 2 end wrenches of 14.

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3.3- How to operate

Important: Before and after every running, turn the tool speed potentiometer to 0.

- Place the Beaker
- Tighten the Beaker with the fixed Beaker and the strap (see photos in annexe).
- Lower the slide and place the turbine to the desired working height.
- Tighten the slide in location with the bakelite plug.
- Adjust the device of positioning of the turbine, with the button 00023271 placed on the left of the column, by aligning the red point on the height value of the used Beaker: For example, for a Beaker with a 200 mm height, place the red point on 200.

Nota: For height of Beaker lower than 170 mm, place the red point on 170 mm. For height of Beaker higher than 270 mm, place the red point on 270 mm.

- Press the button S1.
- Adjust the tool speed with the potentiometer.
- At the end of the mixing, turn the potentiometer on 0.
- Press the button S0.
- Unscrew the slide and lift it in high location.

Important: You must tighten the Beaker with the strap to obtain the running authorisation.

3.4- Emulsifier assembling (Rotor-Stator type)

- Dismantle the spindle the 2 end wrenches of 10 and 14.
- Lift the positioning the stator dog in the hole placed on the motor slack.
- Tighten the main spindle (rotor) with the wrenches.

Warning: The bad setting of the indexing dog can provoke material damages (see photos in annexe).

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4. ADVICE FOR UTILIZATION

4.1- Safety instructions

- The operator will have to be enough qualified to insure the good adjustment of the device of security.
- In the case of projection or splash risks reduce immediately the speed of the tool.
- Never to make turn the spindle of agitation to large void speed.
- Never to make turn the spindle of agitation to low speed (< 300 rpm) during too long.

4.2- Contra-indications

Important: the running of this machine is forbidden in explosive area or flammable.

- In these case, call us.
- During harmful product mix for the man, take all necessary precautions for avoid all contacts. Especially, eliminate all projections or splash risks.

5. MAINTENANCE

Before every operation, switch off the turbotest.

Cleaning:

- Wipe columns regularly with a dry rag
- Clean the rest of the machine with a humid sponge
- After each utilization, clean tools to the hot water, or possibly with a solvent chlorinated, aromatic (toluen, white spirit...) or the acetone.
- Call us for specific applications.
- Never to use abrasive products.

Motor: sweep change

Never to slip the rotor from the stator because of risk of final demagnetization.

- Check once year the sweeps by removing the two plugs placed on the sides motor.
In case if damage:
- Remove the used sweeps.
- Replace the new sweeps.
- Replace the two plugs.

Fuses:

- Remove the plug of the fuse holder, placed on the back side of the control panel, with a screwdriver.
- Change the fuse by respecting imperatively the caliber.
- Replace the plug of the fuse holder.

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6. TROUBLE SHOOTING

The slide is hard to manipulate:

- Check that the bakelite plug placed on the left side of the slide is unscrewed.
- Check the cable.
- Check the cleaning of the guidance columns.

The slide stays in low position:

- Check the cable
- Check the catching points of the cable on the slide and the counter weight.

The tool does not turn:

- Check the location of the red point on the ruler: the read value must be lower or equal than the Beaker height.
- Check the good locking of the Beaker.
- Check the fuse.
- Check the motor sweeps.
- Check the limit switch of the slide

The push button S1 does not switch on any more:

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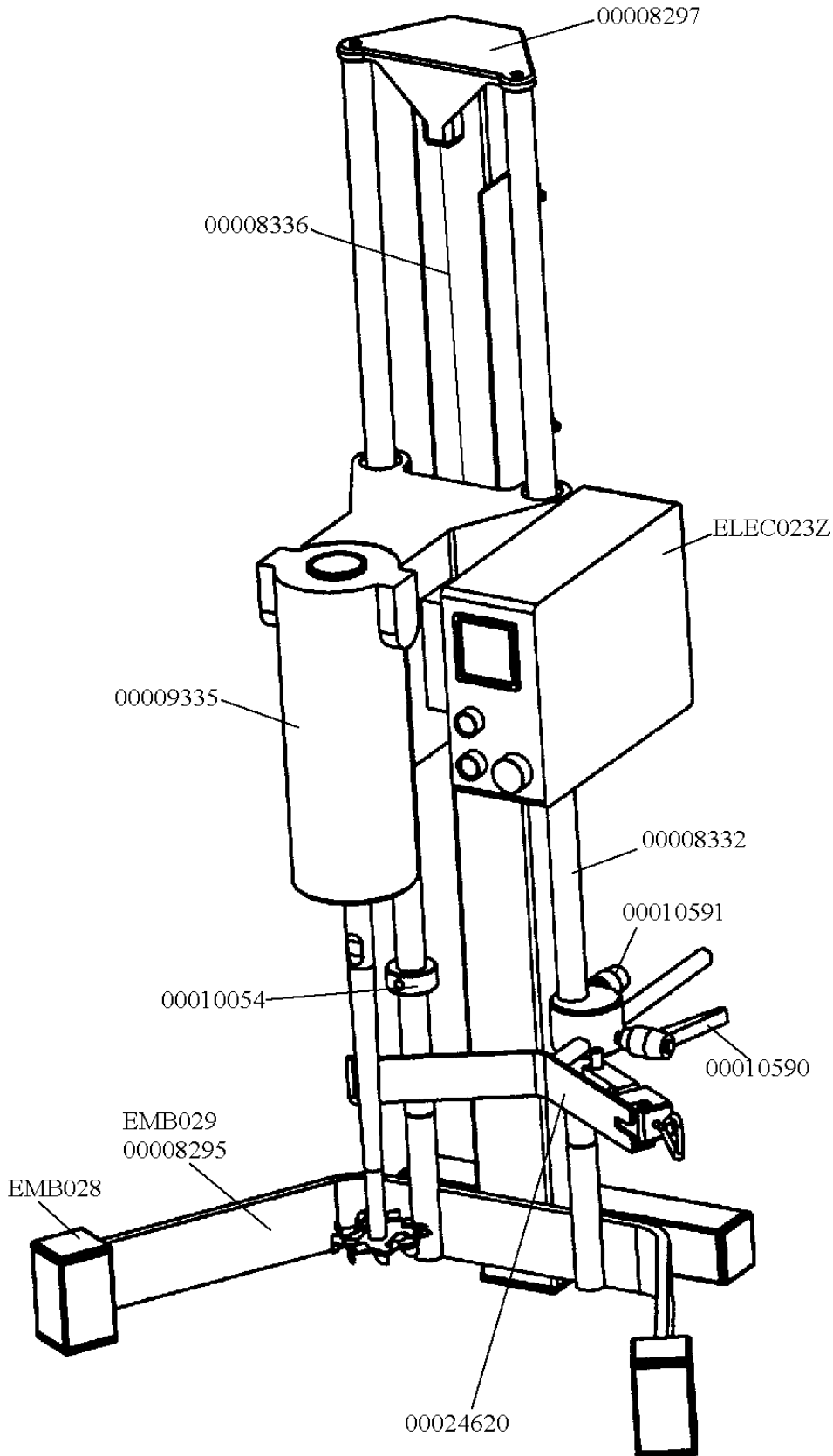
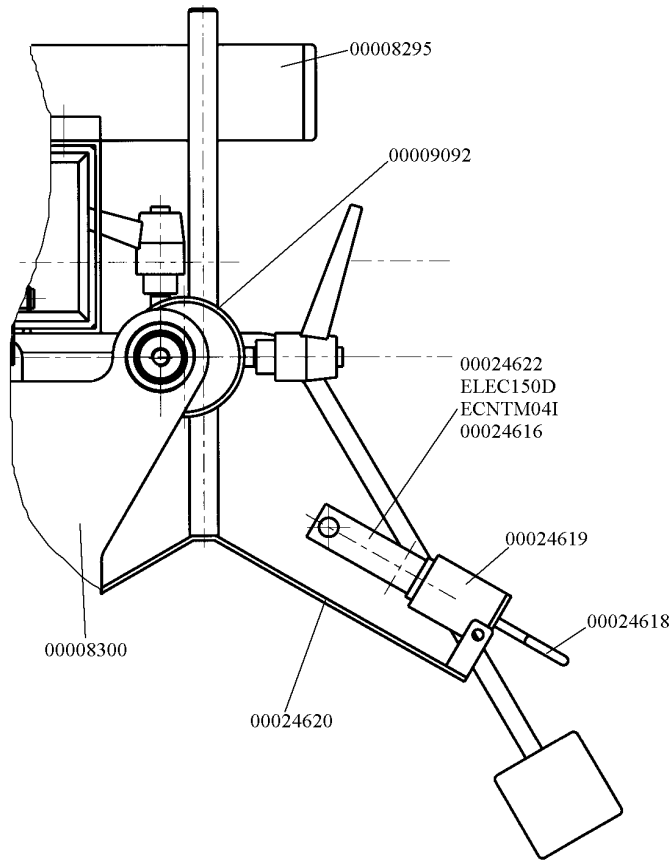
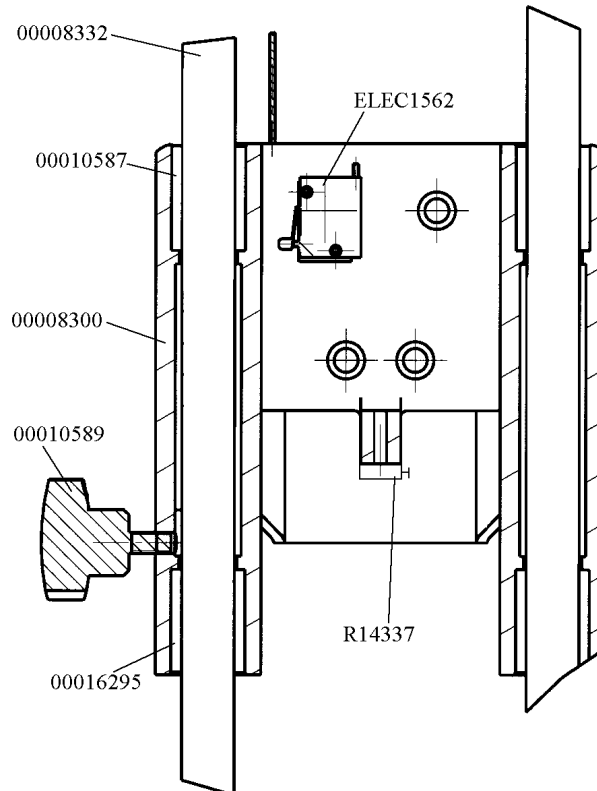
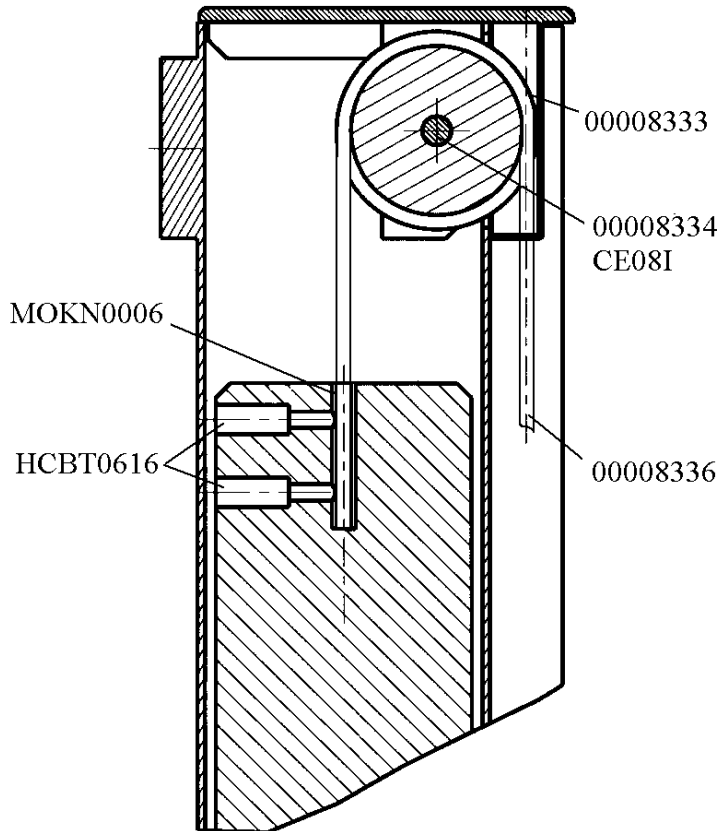
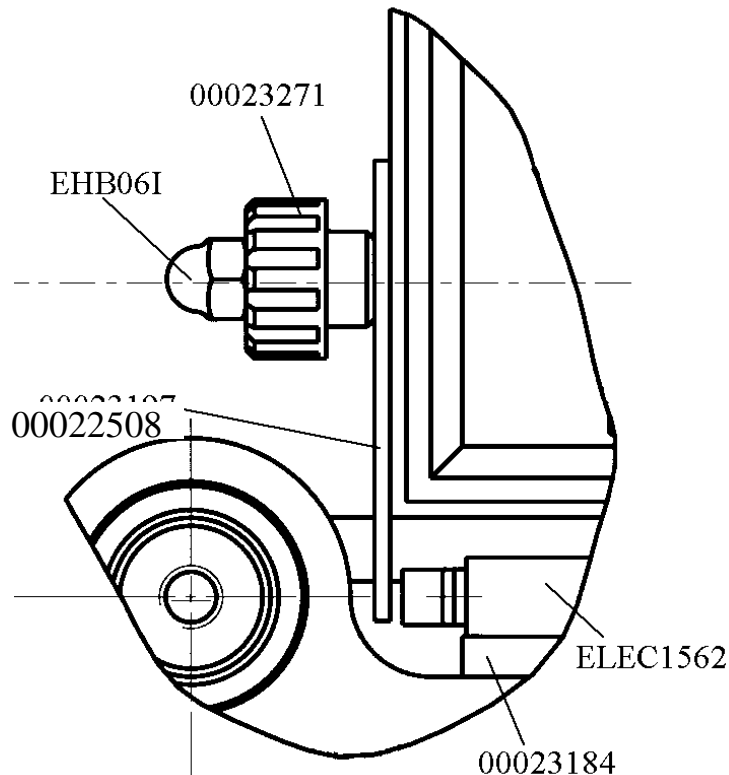
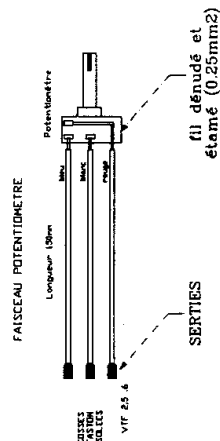
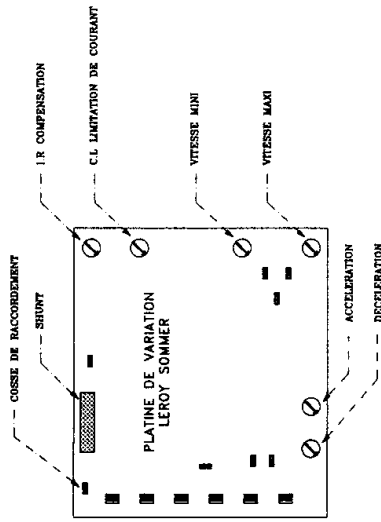
9. ANNEXE

9.1- Conformity

- With the directive 89/392/EEC
- With the working code

9.2- Extract from the conditions of sale

• **Guarantee:**

⇒The normal guarantee period for equipment is 6 months. It can never exceed one year even in special cases which have been explicitly negotiated. In case of equipment which is used day and night, the guarantee is automatically halved.

⇒It is strictly limited to replacements or repairs of parts which no longer work due to faults or construction errors.

⇒It does not apply to replacements or repairs which result from normal use of apparatus or machines, nor to deterioration or accidents caused by negligence, lack of surveillance, servicing or defective installation and use of apparatus.

• **Transports:**

⇒All transport operations, insurance, customs duties, taxes and handling are at the purchaser's charge, as well as the risks and perils resulting from the transportation, and it is the purchaser's duty to verify on arrival of the goods that is well, and if need be to make a claim against the transporters even if the goods have been sent at our charge.

⇒In case of goods being sent by the seller, this is done at the lowest rates possible and the transport is payable on reception of the goods, except if the purchaser has expressly demanded otherwise, and in all cases it is his entire responsibility.

• **Claim:**

⇒Any claim relating to good supplied or their payment are to be treated by the "Tribunal de commerce de Nantes", which is the sole authority no matter what the conditions of sale or where there exists several defendants.

9.3- After sales service

For the maintenance, first call the seller of the machine.

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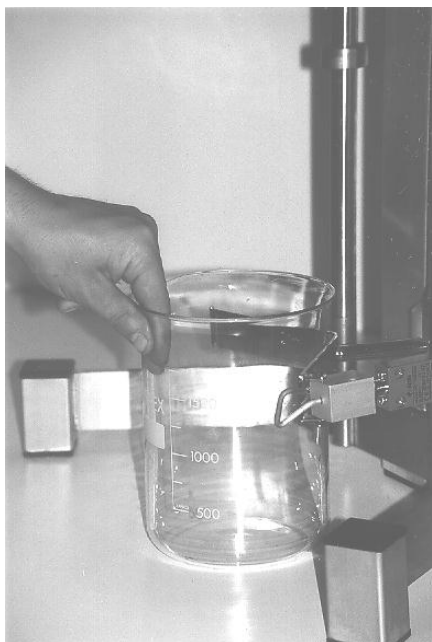
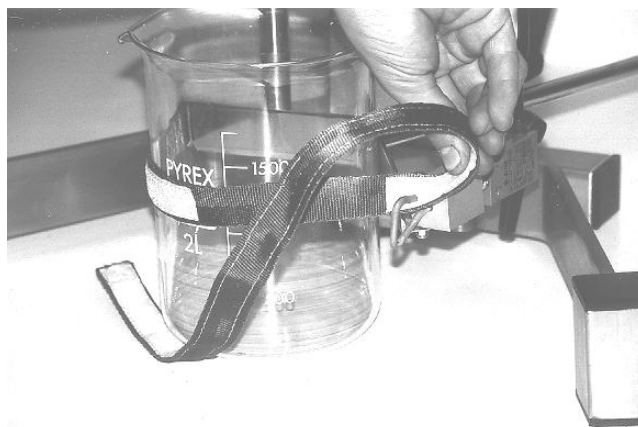
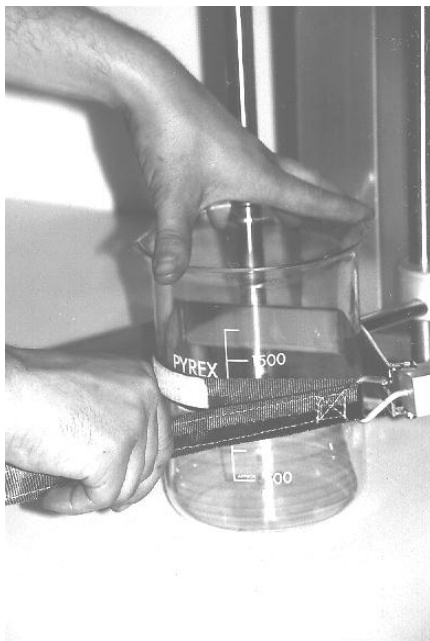
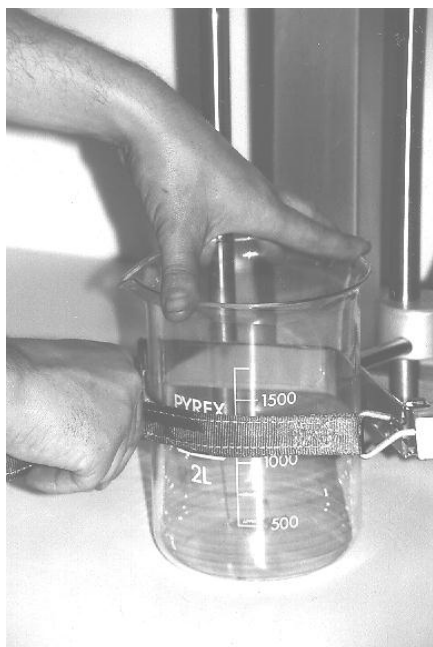
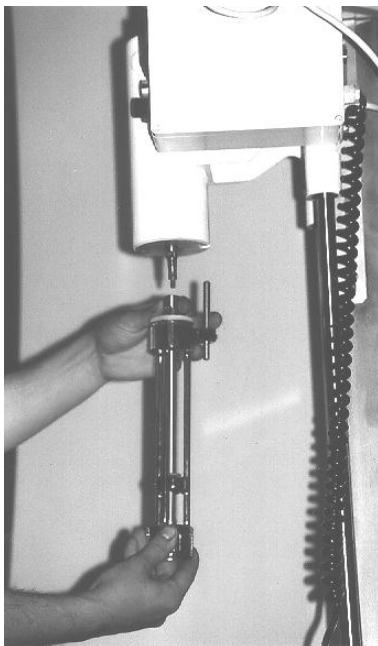
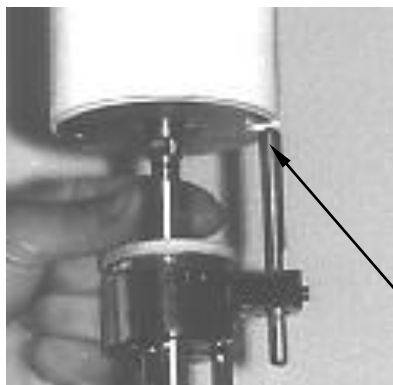
EMULSIFIER ASSEMBLING (indexing dog setting up)

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