English	User Manual Printer P25, P26, P28
Deutsch	Benutzerhandbuch Drucker P25, P26, P28
Español	Manual de usuario Impresora P25, P26 y P28
Français	Guide de l'utilisateur Imprimante P25, P26, P28
简体中文	简明用户手册 打印机 P25, P26, P28
日本語	ユーザマニュアル プリンタ P25, P26, P28

METTLER TOLEDO

This User Manual is a brief instruction that provides information to handle with the first steps of the instrument in a safe and efficient manner. Personnel must have carefully read and understood this manual before performing any tasks.

For full information, always consult the Reference Manual (RM).

► www.mt.com/printer-p-20-RM

Overview printer

1	On/Off Switch	2	Pilot lamp – status indicator
3	Control panel	4	Interface connector
5	Power connection		
1	Ein-/Ausschalter	2	Kontrolllampe – Statusanzeige
3	Bedienelement	4	Schnittstellenanschluss
5	Stromanschluss		
1	Interruptor de encendido/apagado	2	Lámpara de indicación (indicador de estado)
3	Panel de control	4	Conector de interfaz
5	Conexión a la alimentación		
1	Interrupteur Marche/Arrêt	2	Voyant témoin – indicateur d'état
3	Panneau de commande	4	Connecteur d'interface
5	Branchement secteur		
1	双控开关	2	指示灯 – 状态指示灯
3	控制面板	4	接口接头
5	电源连接		
1	オン / オフスイッチ	2	パイロットランプ – ステータスインジケータ
3	制御パネル	4	インターフェースコネクタ
5	電源コネクタ		

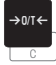
Overview functions

Key	P25	P26	P28
	<ul style="list-style-type: none"> Paper feed Menu option (+ On switch) 	<ul style="list-style-type: none"> Paper feed Scroll up menu option 	<ul style="list-style-type: none"> Paper feed Scroll up menu option
	–	<ul style="list-style-type: none"> Zeroing / taring the balance Abort – exit 	<ul style="list-style-type: none"> Zeroing / taring the balance Abort – exit
	–	<ul style="list-style-type: none"> Printout of date and time Scroll down menu option 	<ul style="list-style-type: none"> Printout of date and time Scroll down menu option
	–	<ul style="list-style-type: none"> Printing of stable weight values Opening the menu – select next menu option Save settings 	<ul style="list-style-type: none"> Printing of stable weight values Opening the menu – select next menu option Save settings
	–	–	<ul style="list-style-type: none"> Totaling application
	–	–	<ul style="list-style-type: none"> Statistics application

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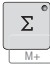
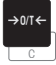
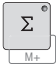
1 Safety Information

Two documents named "User Manual" and "Operating Instructions" are available for this instrument.

- The User Manual shows you how to start using the instrument.
- The Operating Instructions contain a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Operating Instructions. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

1.1 Further applicable documents

This User Manual is a brief instruction that provides information to handle with the first steps of the instrument in a safe and efficient manner. Personnel must have carefully read and understood this manual before performing any tasks.

For full information, always consult the Reference Manual (RM).

► www.mt.com/printer-p-20-RM

1.2 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

WARNING	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
CAUTION	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
NOTICE	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

Warning symbols

Electrical shock

General hazard: read the Operating Instructions for information about the hazards and the resulting measures.

Notice

1.3 Product specific safety notes

Intended use

This instrument is designed to be used in laboratories by trained staff. The instrument is intended for print out protocol purposes.

Any other type of use and operation beyond the limits of technical specifications without written consent from Mettler-Toledo GmbH is considered as not intended.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

METTLER TOLEDO assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. METTLER TOLEDO assumes that the instrument owner provides the necessary protective gear.

Safety notes

WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the approved METTLER TOLEDO power supply cable and AC/DC adapter with a current-limited SELV output.
- 2 Connect the power cable to a grounded power outlet, ensure correct polarity.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and power plug for damage and replace damaged cables and power plugs.

NOTICE

Damage to the instrument due to the use of unsuitable parts

Using unsuitable parts with the instrument can damage the instrument or cause it to malfunction.

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

2 Installation and Putting into Operation

2.1 Scope of delivery

- AC/DC adapter
- Power cable (country specific)
- User Manual
- Declaration of conformity

2.2 Connecting the printer to the power supply

WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the approved METTLER TOLEDO power supply cable and AC/DC adapter with a current-limited SELV output.
- 2 Connect the power cable to a grounded power outlet, ensure correct polarity.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and power plug for damage and replace damaged cables and power plugs.

NOTICE

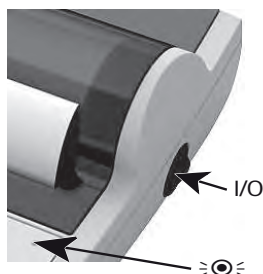
Damage to the AC/DC adapter due to overheating

If the AC/DC adapter is covered or in a container, it is not sufficiently cooled and will overheat.

- 1 Do not cover the AC/DC adapter.
- 2 Do not put the AC/DC adapter in a container.

The printer is supplied with a universal AC/DC adapter or an AC/DC adapter with a country-specific power cable.

- Install the cables so that they cannot be damaged or interfere with operation.
 - Insert the power cable in a grounded power outlet that is easily accessible.
- 1 Connect the AC/DC adapter to the connection socket on the back of the printer.
 - 2 Lock the connector with the screw nut.
 - 3 Connect the 3-pin grounded power cable to the socket for the AC/DC adapter.
 - 4 Connect the printer and the balance with the interface cable.
 - 5 Switch the balance on.
 - 6 Connect the printer to the power supply.
 - 7 Switch on the printer with the **I/O** switch.

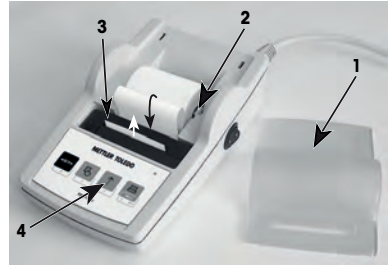
 **Note**

RS232 interface only; the pilot lamp will flash until an instrument is connected. If the lamp is still flashing check the chapter "Troubleshooting" in the Reference Manual.

2.3 Inserting paper

- 1 Remove paper cover (pull on the back) (1).
- 2 Insert the paper rotary axis through the core of the paper roll (2).
- 3 Feed paper through the slot in the printing unit in the direction of the arrow (3).
- 4 Press and hold until enough paper has been fed (4).
- 5 Lead paper through the paper cover slot.
- 6 Place the paper cover (1) on.

If, after the first printouts the connection settings (Bauderate, Bit, Handshake) of the balance change, the printer must be restarted. Switch off / on to start the auto bauderate detection again. Auto bauderate can be switched off.

3 Keystroke Functions

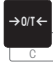
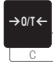
Printer P25

Operation level	 Press briefly	 Press and hold down
	one line feed	continuous feed (until release)
 +  On	Printer test with setting printout	open menu
Menu level	 Press briefly	 Press and hold down
	next menu item	store

Printer P26

Operation level	 Press briefly	 Press and hold down		
	zero / tare the balance	–		
	print Date and Time	–		
	one line feed	continuous feed (until release)		
	print	open menu		
Menu level	 Press briefly	 Press and hold down		
 exit	 down	 up	 next	 store

Printer P28

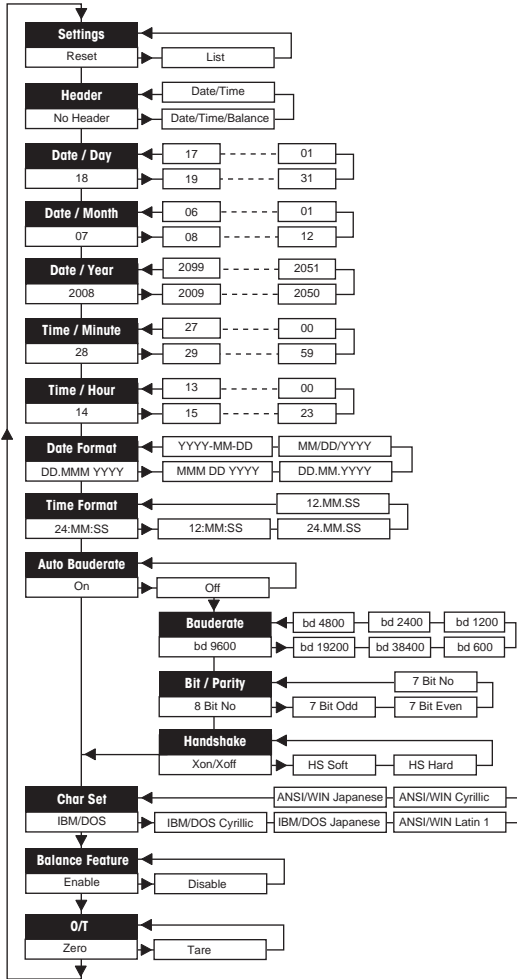
Operation level	 Press briefly	 Press and hold down
	zero / tare the balance	–
	print Date and Time	–
	one line feed	continuous feed (until release)
	print	open menu
Menu level	 Press briefly	 Press and hold down
 exit  down  up  next	 store	
Function level	 Press briefly	 Press and hold down
	Start totaling application	Conclude totaling application
	Start statistics application	Conclude statistics application

4 Menu

4.1 Printer P25

Press briefly	Press and hold down	Printout
<div style="border: 1px solid black; padding: 2px; text-align: center;">Settings</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Reset</div>		Settings Reset
<div style="border: 1px solid black; padding: 2px; text-align: center;">Reset the printer</div>		<pre> ----- Reset done ----- Current settings: Baudrate: 9600 Bit/Parity: 8 Bit Non Handshake: Off Char Set: IBM/DOS </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">RS232 interface only</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Auto Bauderate</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">On</div>	Store	<pre> Auto Bauderate: On -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">Auto Bauderate</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Off</div>	Store	<pre> Auto Bauderate: Off -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">Char Set</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">IBM / DOS</div>	Store	<pre> IBM/DOS -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">IBM / DOS</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Cyrillic</div>	Store	<pre> IBM/DOS Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">IBM / DOS</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Japanese</div>	Store	<pre> IBM/DOS Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Latin 1</div>	Store	<pre> ANSI/WIN Latin 1 -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Cyrillic</div>	Store	<pre> ANSI/WIN Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Japanese</div>	Store	<pre> ANSI/WIN Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">Balance Feature</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Enable</div>	Store	<pre> Balance Feature: Enable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">Balance Feature</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Disable</div>	Store	<pre> Balance Feature: Disable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 2px; text-align: center;">Menu Option</div> <div style="border: 1px solid black; padding: 2px; text-align: center;">Factory Setting</div>		

4.2 Printer P26 / P28

Examples if balance display is used:

Settings	Reset	<i>rESEt</i>
	List	<i>LlSt</i>
Header	*No Header	<i>no Hdr</i>
	Date/Time	<i>d t Hdr</i>
	Date/Time/Balance	<i>dt b Hdr</i>
Date Day**	18	<i>dRY 18</i>
Date Month**	07	<i>Month 07</i>
Date Year**	2008	<i>YEAR-2008</i>
Time Minute**	28	<i>Min. 28</i>
Time Hour**	14	<i>Hour 14</i>
Date Format***	DD.MMM.YYYY	<i>ddMMn Y</i>
	MM/DD/YYYY	<i>MM dd Y</i>
	DD.MM.YYYY	<i>ddMMY</i>
	MM/DD/YYYY	<i>MM dd Y</i>
	YYYY-MM-DD	<i>Y-MM-dd</i>
Time Format***	24:MM:SS	<i>24:MM:SS</i>
	12:MM:SS	<i>12:MM:SS</i>
	24.MM.SS	<i>24MMSS</i>
	12.MM.SS	<i>12MMSS</i>
Baud rate	*9600	<i>bd9600</i>
	19200	<i>bd.19200</i>
	38400	<i>bd.38400</i>
	600	<i>bd600</i>
	1200	<i>bd.1200</i>
	2400	<i>bd.2400</i>
	4800	<i>bd.4800</i>
Bit/Parity	*8 Bit No	<i>8b-no</i>
	7 Bit Odd	<i>7b-odd</i>
	7 Bit Even	<i>7b-e</i>
	7 Bit No	<i>7b-no</i>
Handshake	*Xon/Xoff	<i>HS off</i>
	HS Soft	<i>HS Soft</i>
	HS Hard	<i>HS Hard</i>
Char Set	*IBM/DOS	<i>dOS</i>
	IBM/DOS Cyrillic	<i>dOS Cyr</i>
	IBM/DOS Japanese	<i>dOS JRP</i>
	Ansi/Win Latin 1	<i>11 in LRE</i>
	Ansi/Win Cyrillic	<i>11 in Cyr</i>
	Ansi/Win Japanese	<i>11 in JRP</i>
Tare/Zero	*Zero	<i>ZErO</i>
	Tare	<i>tAR-E</i>

Menu Option
Factory Setting

* factory settings

** instruments with built-in clock: not applicable, built-in clock has priority

*** instruments with built-in clock: format of printer has priority

Example for setting the day (P26 / P28 only)

Open the menu

- Hold down the  key until **Reset** appears.
- Select the menu option **Day** with the  key.
- Select the day, e.g. day 21.
Scroll up with the \uparrow key.
Scroll down with the \downarrow key.

Save the settings

- Hold down the  key until **Stored** appears.

Statistics application (P28 only)

Sample weight must be in the range of 70% - 130% of the current average value.
Minimum dynamic weight changed must be 100 d.

- 1 Press  to start the Statistics application (LED on).
- 2 Add a new sample and press **M+**.
- 3 Press  to print subtotal, if needed.
- 4 Press and hold  to print and end the Statistics application (LED off).

Note

- Unit change during the application is not permitted.
- Maximal 999 sample.

Totaling application (P28 only)

Minimum weights + 1 d.

- 1 Press  to start the Totaling application (LED on).
- 2 Add a new sample and press **M+**.
- 3 Press  to print subtotal, if needed.
- 4 Press and hold  to print and end the Totaling application (LED off).

Note

- Unit change during the application is not permitted.
- Maximal 999 sample.

5 Maintenance

WARNING

Death or serious injury due to electric shock

Contact with parts carrying a live current can lead to injury and death.

- 1 Disconnect the printer from the power supply prior to cleaning and maintenance.
- 2 Prevent liquid from entering the printer or AC/DC adapter.

5.1 Replacing the paper roll

The insertion of the new paper roll is described in chapter [Inserting paper ▶ Page 6]. The order number for the paper roll can be found in chapter [Consumables ▶ Page 14].

Note

The paper rotary axis must always be transferred from the old paper roll to the new one.

5.2 Replacing the ribbon

The order number for the ribbon cartridge can be found in chapter [Consumables ▶ Page 14].

- 1 Remove the paper cover (1).
- 2 Pull paper out of printing unit (2).

- 3 Remove ribbon cartridge, by pressing in arrow direction (3).
- 4 Insert the new ribbon cartridge. If necessary, take up slack with tensioning wheel (4).

Insertion of a new paper roll is described in chapter [Inserting paper ▶ Page 6].

5.3 Cleaning the housing

Since the printer housing is made of top grade, resistant materials, all commercially available mild cleaning agents may be used.

NOTICE

Damage to printer due to inappropriate cleaning methods

The printer housing is made from high quality, resistant materials and can be damaged by certain cleaning agents, solvents or abrasives. Any liquid that enters the housing may damage the printer.

- 1 Use water and a mild detergent to clean the printer.
- 2 Wipe off any spills immediately.
- 3 Prevent liquid from entering the interior of the printer.

Contact a METTLER TOLEDO representative to find about the service options available – regular maintenance by an authorized service engineer will ensure consistent weighing accuracy over the long term and extend the service life of the instrument.

6 Technical Data

6.1 General technical data

Power supply

AC/DC adapter:	Primary: 100 – 240 V, 50/60 Hz Secondary: 12 V DC, 2.5 A (with electronic overload protection)
Cable for AC/DC adapter:	3-core, with country-specific plug
Polarity:	 with a current limited SELV (Safety Extra Low Voltage) output
Power line voltage:	12 V DC
Power consumption:	Max. 2.5 A

Protection and standards

Overvoltage category:	II
Degree of pollution:	2
Standards for safety and EMC:	See Declaration of Conformity
Range of application:	For use in closed interior rooms only

Environmental conditions

Height above mean sea level:	Up to 4000 m
Ambient temperature:	0 to 40 °C
Storage condition:	-25 to 60 °C
Relative air humidity:	Max. 80% up to 31 °C, linearly decreasing to 50% at 40 °C, noncondensing

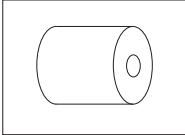
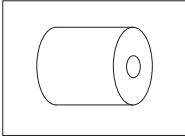
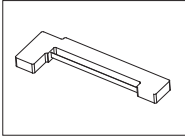
Printer features

Print technology:	Dot matrix printer 5 × 7, 24-character line length
Print speed:	2.3 lines per second
Ribbon cartridge:	Exchangeable, black
Paper roll:	Standard paper 57.5 mm × ø 50 to 60 mm, integrated in housing, commercial size
Print quality:	Light resistant and thermally stable printing (GLP, GMP, ISO 9001)
Interfaces:	RS-P25, RS-P26, RS-P28 USB-P25
Compatibility limits:	RS232C USB <ul style="list-style-type: none"> Function Date / Time: not applicable on balances with built-in clock (built-in clock has priority) Function Zeroing / Taring: not applicable on all models of balances

Dimension / weight

Printer dimensions (W × D × H):	120 × 203 × 73 mm
Packaging dimensions (W × D × H):	255 × 205 × 150 mm
Net weight:	740 g (paper roll incl.)

7 Consumables

Description	Part No.	
 A line drawing of a cylindrical paper roll, shown from a perspective view. The roll is oriented horizontally, with a circular end face on the right side.	Paper roll (length: 20 m), set of 5 pcs	00072456
 A line drawing of a cylindrical paper roll, shown from a perspective view. The roll is oriented horizontally, with a circular end face on the right side.	Paper roll (length: 13 m), self-adhesive, set of 3 pcs	11600388
 A line drawing of a ribbon cartridge, shown from a perspective view. It is a long, thin, rectangular component with a small protrusion on the left side and a notch on the right side.	Ribbon cartridge, black, set of 2 pcs	00065975

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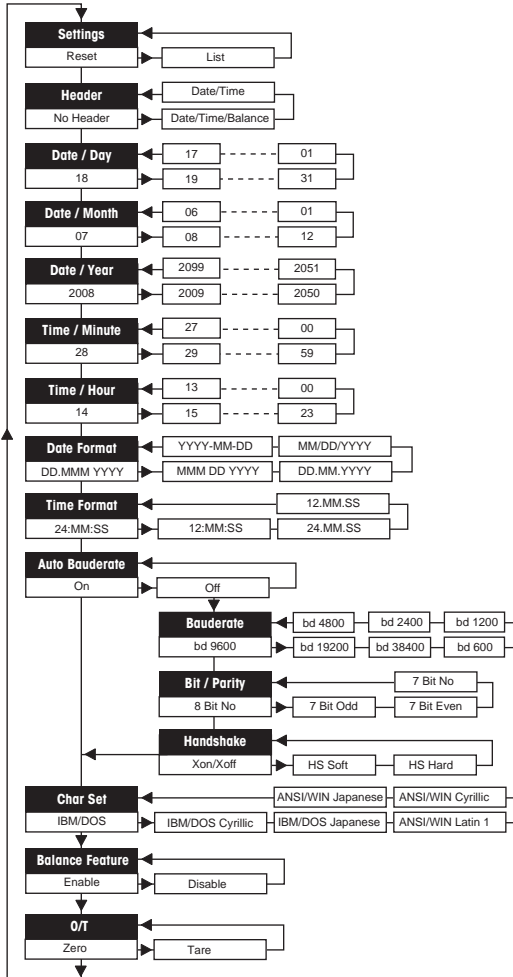
4 Menü

4.1 Drucker P25

Kurz drücken	Gedrückt halten	Ausdruck
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Settings Reset </div>		Settings Reset
	Drucker zurücksetzen	<pre> ----- Reset done ----- Current settings: Baudrate: 9600 Bit/Parity: 8 Bit Non Handshake: Off Char Set: IBM/DOS </pre>
<div style="border: 1px solid gray; padding: 5px; text-align: center;"> Nur RS232-Schnittstelle <div style="border: 1px solid black; padding: 5px; text-align: center;"> Auto Bauderate On </div> </div>	Speichern	<pre> Auto Bauderate: On -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid gray; padding: 5px; text-align: center;"> <div style="border: 1px solid black; padding: 5px; text-align: center;"> Auto Bauderate Off </div> </div>	Speichern	<pre> Auto Bauderate: Off -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Char Set IBM / DOS </div>	Speichern	<pre> IBM/DOS -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> IBM / DOS Cyrillic </div>	Speichern	<pre> IBM/DOS Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> IBM / DOS Japanese </div>	Speichern	<pre> IBM/DOS Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Latin 1 </div>	Speichern	<pre> ANSI/WIN Latin 1 -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Cyrillic </div>	Speichern	<pre> ANSI/WIN Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Japanese </div>	Speichern	<pre> ANSI/WIN Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Balance Feature Enable </div>	Speichern	<pre> Balance Feature: Enable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Balance Feature Disable </div>	Speichern	<pre> Balance Feature: Disable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Menüoption Werkseinstellung </div>		

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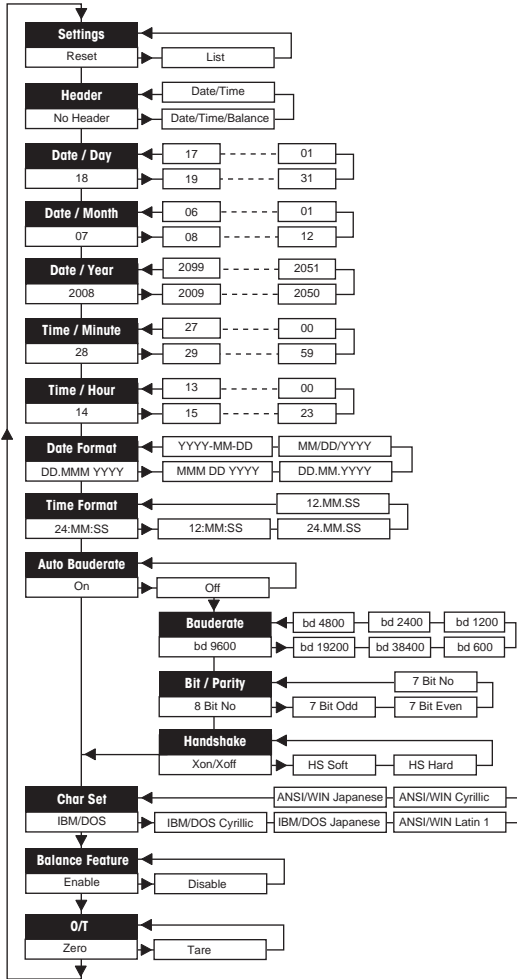
4 Menú

4.1 Impresora P25

Pulsación breve	Pulsación prolongada	Impresión
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Settings Reset </div>		Settings Reset
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Solo para la interfaz RS232 Auto Bauderate On </div>	Reiniciar la impresora	<pre> ----- Reset done ----- Current settings: Baudrate: 9600 Bit/Parity: 8 Bit Non Handshake: Off Char Set: IBM/DOS </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Solo para la interfaz RS232 Auto Bauderate Off </div>	Guardar	<pre> Auto Bauderate: On -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Auto Bauderate Off </div>	Guardar	<pre> Auto Bauderate: Off -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Char Set IBM / DOS </div>	Guardar	<pre> IBM/DOS -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> IBM / DOS Cyrillic </div>	Guardar	<pre> IBM/DOS Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> IBM / DOS Japanese </div>	Guardar	<pre> IBM/DOS Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Latin 1 </div>	Guardar	<pre> ANSI/WIN Latin 1 -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Cyrillic </div>	Guardar	<pre> ANSI/WIN Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> ANSI / WIN Japanese </div>	Guardar	<pre> ANSI/WIN Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Balance Feature Enable </div>	Guardar	<pre> Balance Feature: Enable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> Balance Feature Disable </div>	Guardar	<pre> Balance Feature: Disable -----STORED----- ... "current settings" </pre>

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24.MM.SS 24-MM-SS
12.MM.SS 12-MM-SS
24.MM.SS 24MMSS
12.MM.SS 12MMSS

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*9600 bd9600
19200 bd.19200
38400 bd.38400
600 bd600
1200 bd.1200
2400 bd.2400
4800 bd.4800

Bit/paridad

*8 Bit No 8b-no
7 Bit Odd 7b-odd
7 Bit Even 7b-e
7 Bit No 7b-no

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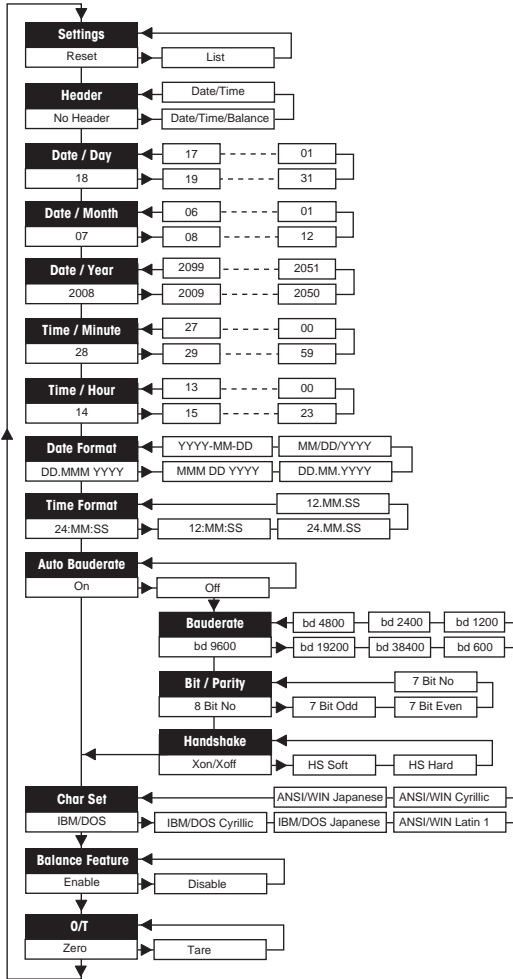
4 Menu

4.1 Imprimante P25

Appuyez brièvement	Appuyez et maintenez la touche enfoncée.	Ticket d'impression
<div style="border: 1px solid black; padding: 5px; text-align: center;">Settings</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Reset</div>		Settings Reset
	Réinitialiser l'imprimante	<pre> ----- Reset done ----- Current settings: Baudrate: 9600 Bit/Parity: 8 Bit Non Handshake: Off Char Set: IBM/DOS </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Interface RS232 uniquement</div>		
<div style="border: 1px solid black; padding: 5px; text-align: center;">Auto Bauderate</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">On</div>	Enregistrer	<pre> Auto Bauderate: On -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Auto Bauderate</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Off</div>	Enregistrer	<pre> Auto Bauderate: Off -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Char Set</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">IBM / DOS</div>	Enregistrer	<pre> IBM/DOS -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">IBM / DOS</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Cyrillic</div>	Enregistrer	<pre> IBM/DOS Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">IBM / DOS</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Japanese</div>	Enregistrer	<pre> IBM/DOS Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Latin 1</div>	Enregistrer	<pre> ANSI/WIN Latin 1 -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Cyrillic</div>	Enregistrer	<pre> ANSI/WIN Cyrillic -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Japanese</div>	Enregistrer	<pre> ANSI/WIN Japanese -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Balance Feature</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Enable</div>	Enregistrer	<pre> Balance Feature: Enable -----STORED----- ... "current settings" </pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Balance Feature</div> <div style="border: 1px solid black; padding: 5px; text-align: center;">Disable</div>	Enregistrer	<pre> Balance Feature: Disable -----STORED----- ... "current settings" </pre>

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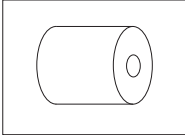
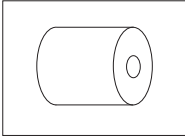
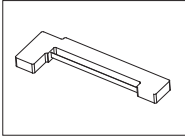
7 Consommables

Description	Référence	
 A line drawing of a roll of paper, shown from a perspective view. It is a cylinder with a smaller circle on the right end representing the core.	Rouleau de papier (longueur : 20 m), 5 unités	00072456
 A line drawing of a roll of adhesive paper, shown from a perspective view. It is a cylinder with a smaller circle on the right end representing the core.	Rouleau de papier adhésif (longueur : 13 m), 3 unités	11600388
 A line drawing of a black ribbon cartridge. It is a long, thin rectangular component with a small notch on the left side and a slightly raised section on the right side.	Cartouche de ruban, noir, 2 unités	00065975

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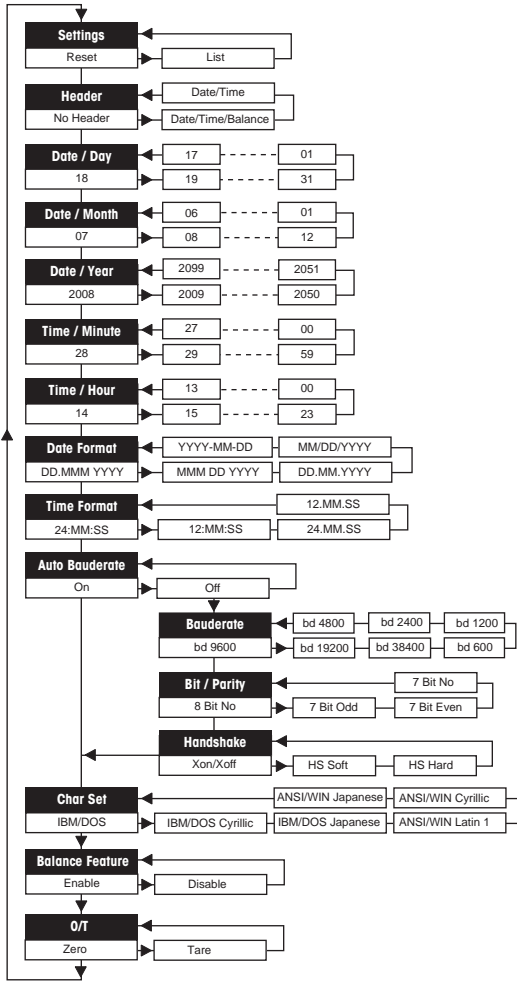
☞ 短按	☞ 向下按住	打印输出
<div style="border: 1px solid black; padding: 5px; text-align: center;">Settings Reset</div>		Settings Reset
	复位打印机	<pre>----- Reset done ----- Current settings: Baudrate: 9600 Bit/Parity: 8 Bit Non Handshake: Off Char Set: IBM/DOS</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;"> 仅限 RS232接 <input type="checkbox"/> </div>		
<div style="border: 1px solid black; padding: 5px; text-align: center;">Auto Bauderate On</div>	存储	<pre>Auto Bauderate: On -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Auto Bauderate Off</div>	存储	<pre>Auto Bauderate: Off -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Char Set IBM / DOS</div>	存储	<pre>IBM/DOS -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">IBM / DOS Cyrillic</div>	存储	<pre>IBM/DOS Cyrillic -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">IBM / DOS Japanese</div>	存储	<pre>IBM/DOS Japanese -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN Latin 1</div>	存储	<pre>ANSI/WIN Latin 1 -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN Cyrillic</div>	存储	<pre>ANSI/WIN Cyrillic -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">ANSI / WIN Japanese</div>	存储	<pre>ANSI/WIN Japanese -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Balance Feature Enable</div>	存储	<pre>Balance Feature: Enable -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">Balance Feature Disable</div>	存储	<pre>Balance Feature: Disable -----STORED----- ... "current settings"</pre>
<div style="border: 1px solid black; padding: 5px; text-align: center;">菜单选项 出厂设置</div>		

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© Mettler-Toledo GmbH 12/2018